


PHARMAGARDES NU-PR797
POSITIVE PRESSURE
COMPOUNDING ASEPTIC ISOLATOR

SLEEVE CHANGING PROCEDURE


1

- [A] Remove Band Clamp
- [B] Remove Sleeve Protector
- [C] Remove Sleeve & Glove Assembly From Isolator

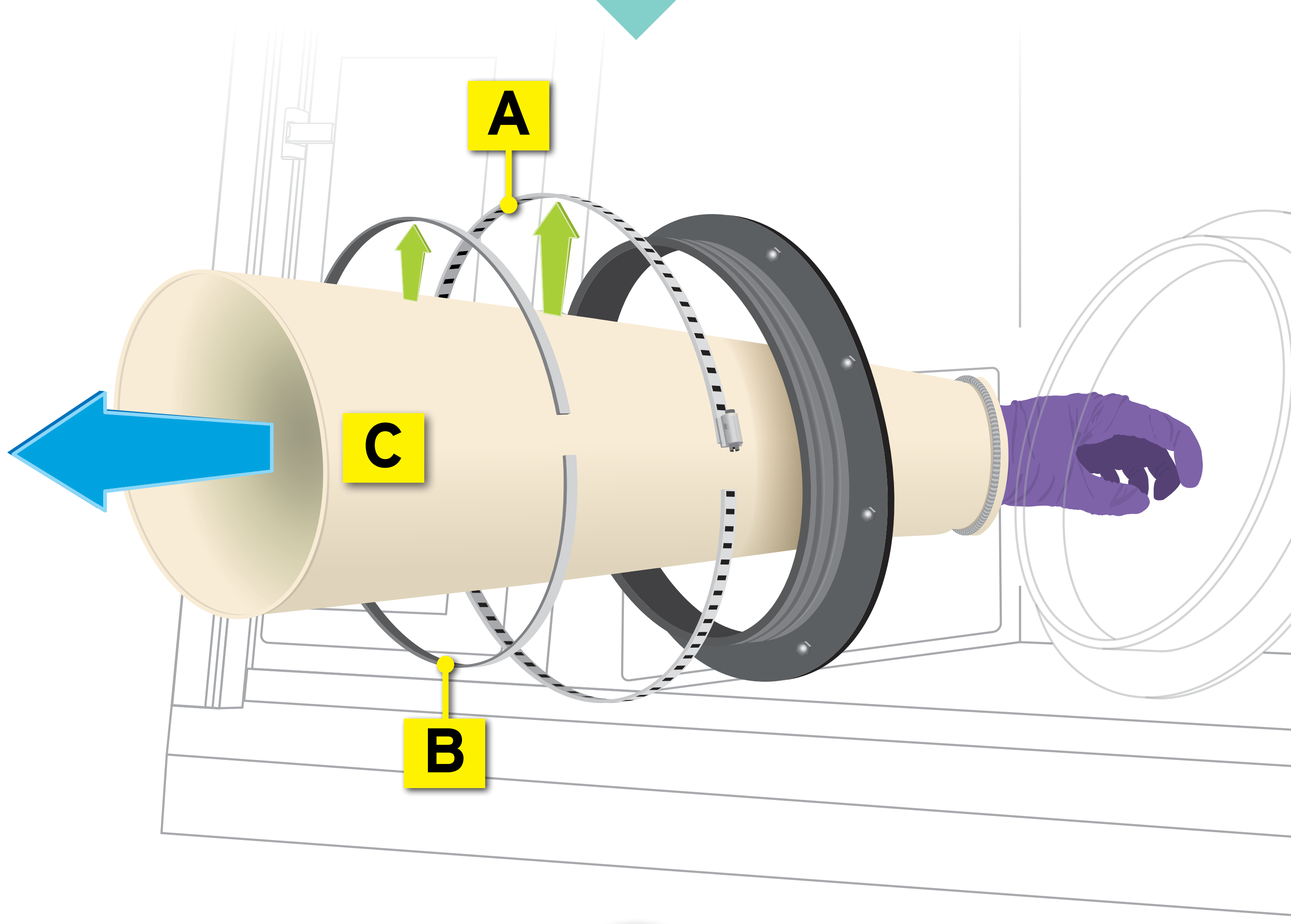


2

- [D] Insert New Sleeve & Glove Assembly Into Isolator
- [E] Roll End of Sleeve Outward and Seat Ring in Second Groove

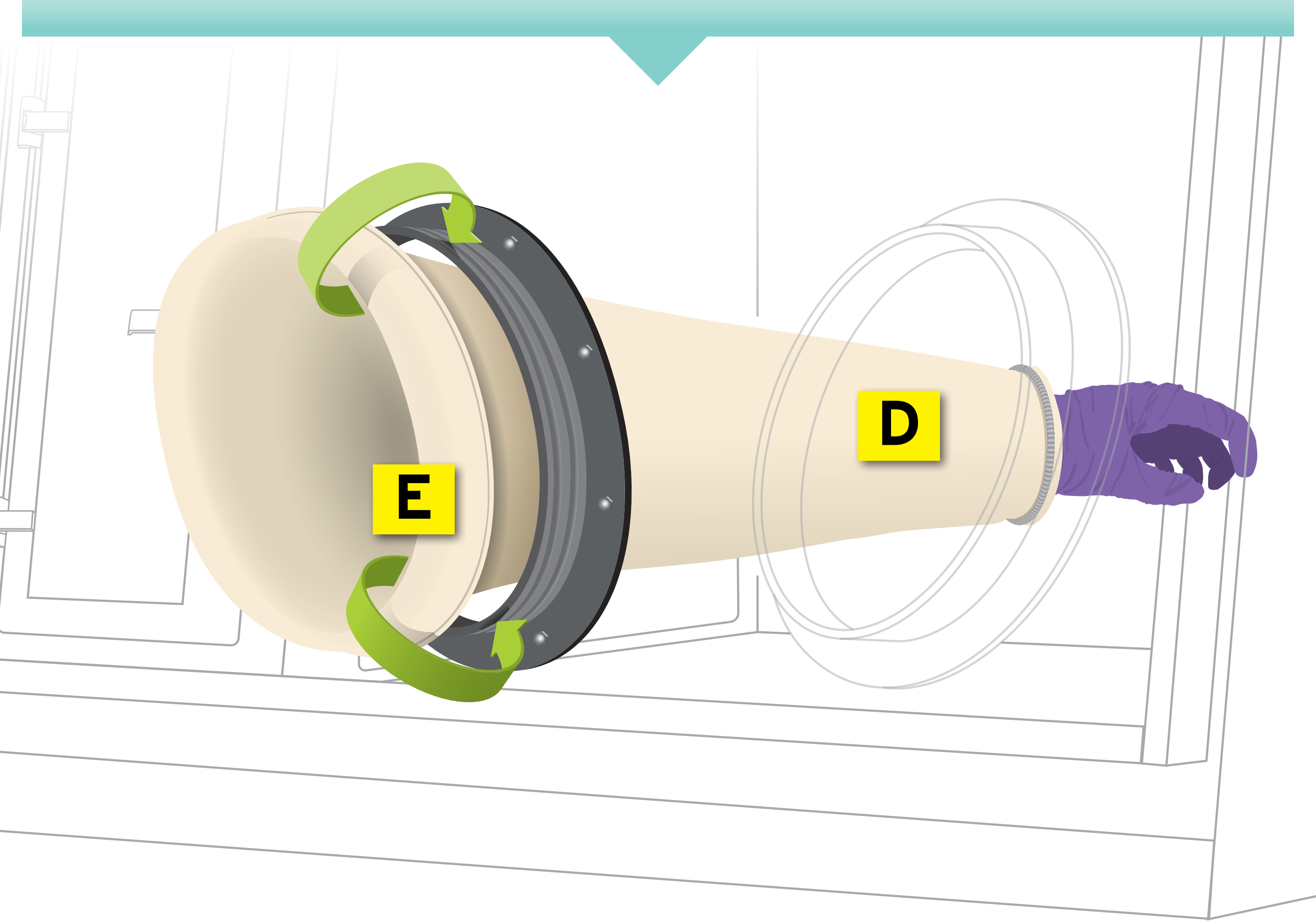

3

- [F] Secure Band Clamp & Sleeve Protector Around New Sleeve

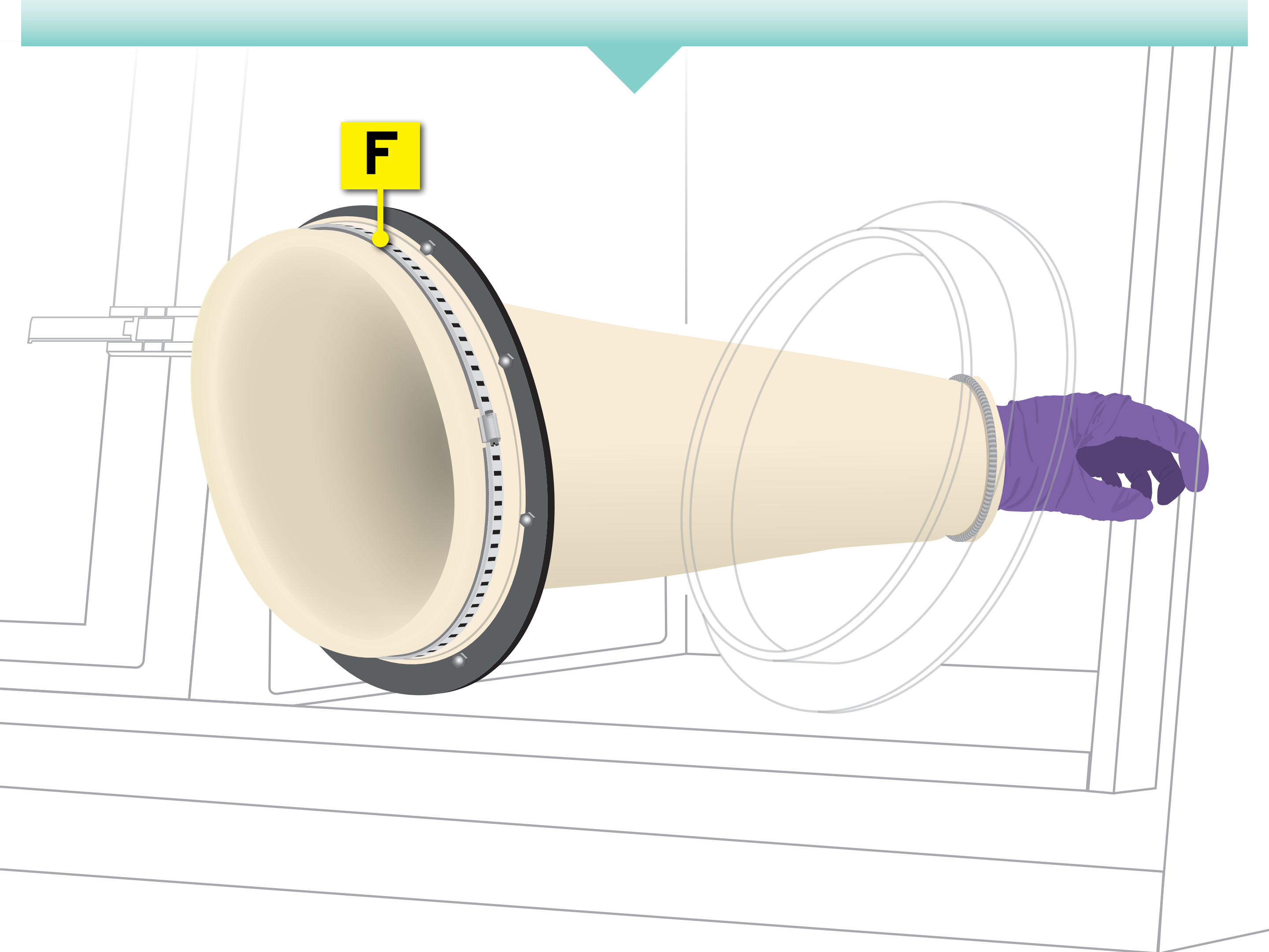

PHARMAGARDES



NuAire, Inc. | nuaire.com