
Good technique when working within an Animal Transfer Station helps protect the animal and limit the spread of allergens.

KNOW THE AIR FLOW AND LOCATION OF AIR CURTAINS

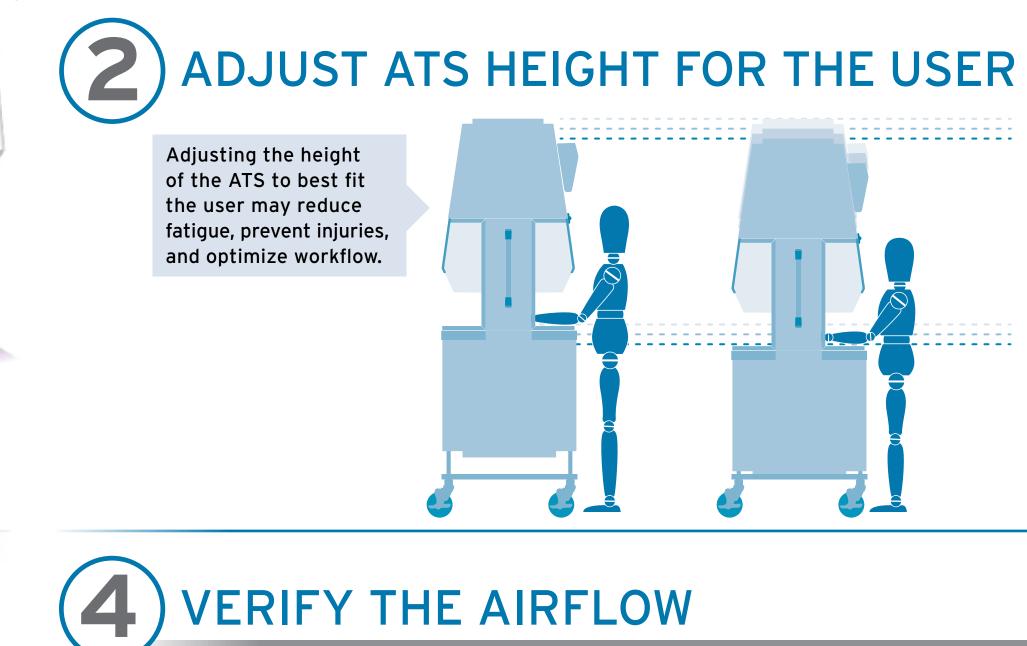

FOR WORKING SAFELY IN YOUR

ANIMAL TRANSFER STATION (ATS)

3) DISINFECT THE WORKSURFACE

Follow disinfecting SOP before

turning on the blower.

Use a disinfectant compatible with surface materials & follow manufacturer's directions. Check the control panel to ensure proper airflow through filters.

Low airflow may indicate that filters need to be cleaned or changed.

NEVER BLOCK THE GRILL ONLY OPEN ONE CAGE WITH ANIMALS AT A TIME Covering the grill can disrupt the air curtain. Only one cage with animals should be open at a time to prevent cross contamination. Follow your facility SOP for handling soiled cages. **POSITION SUPPLIES PREPARE THE CLEAN CAGE FIRST AND WORK FROM CLEAN TO SOILED** IN OR NEAR THE ATS **ON THE CLEAN SIDE** When possible, maintain a clean side and a soiled side in the ATS. Clean cages may be positioned on a cart adjacent to the BSC. Follow SOPs for soiled cages.

