

**TECHNICAL BULLETIN  
ACCESSORY INFORMATION**



2100 Fernbrook Lane  
Plymouth, MN 55447-4722  
U.S.A.

**NU-797 DATA/POWER CORD  
PASS-THRU OPTION (797M383)  
INSTALLATION INSTRUCTIONS**

Phone: 763.553.1270  
Fax: 763.553.0459  
Toll Free: 1.800.328.3352

The NuAire NU-797 data/power cord pass-thru option is designed to allow customers to pass any data or power cord cable through the wall of the Isolator. The data/power cord pass-thru option utilizes two layers of silicone gasket allowing the larger connector cord to be pushed through the Isolator wall. Then the cord is permanently sealed in place with clear silicone RTV. Replacement gaskets and sealers may be purchased as a retrofit kit (RF436) if cable replacement is necessary.



Pass cord through port and position as will be permanently used. NOTE, once cord is sealed into position, no cord movement can take place.



Fill area between inside and outside gaskets with clear silicone RTV provided. Smooth out any excess silicone around the gasket and cord. Allow to set for at least 24 hrs.