



| Parts List | | | | | | |
|------------|-----|---------------|-------|-------|-------|---|
| ITEM | QTY | PART NUMBER | 3 FT. | 4 FT. | 6 FT. | DESCRIPTION |
| 1 | 2 | A-425-7373 | | | | BRACE LATERAL, BSC |
| 2 | 8 | A-500-395 | | | | BUSHING, BASE STAND LEG |
| 3 | 2 | B-400-9562 | | | | WELDMENT, END PANEL, BASESTAND (28") |
| 4 | 4 | B-500-390 | | | | LEG EXTENSION, BASE STAND |
| 5 | 1 | B-500-391-TAB | -01 | -02 | -04 | WELDMENT, CROSS MEMBER PANEL 3 FT. |
| 6 | *4 | B-800-1579 | | | | LEG EXTENSION, BASE STAND EXTENDED HEIGHT |
| 7 | 1 | B-813-1510-01 | | | | BRACKET, 813 BASE STAND, FRONT LEFT |
| 8 | 1 | B-813-1510-02 | | | | BRACKET, 813 BASE STAND FRONT RIGHT |
| 9 | 1 | B-813-1511-01 | | | | BRACKET, 813 BASE STAND, REAR LEFT |
| 10 | 1 | B-813-1511-02 | | | | BRACKET, 813 BASE STAND, REAR RIGHT |
| 11 | 4 | X-980443 | | | | LEG LEVELER FOR 2" SQ TUBING |
| 12 | 8 | X-980493 | | | | ACORN NUT, 1/4-20 |
| 13 | 8 | X-980494-03 | | | | NUT, ACORN 3/8-16 18-8 SST |
| 14 | 8 | X-980495 | | | | 1/4-20 NUTSERT |
| 15 | 12 | X-980604-02 | | | | BOLT, G2, 1/4-20 X .75 LG |
| 16 | 4 | X-980609-06 | | | | SCR PPH 8-32 X 1-1/4 SST |
| 17 | 8 | X-980659-03 | | | | SCREW, HHCS, 1/4-20 X 1/2 18-8 SST |
| 18 | 8 | X-980781-07 | | | | 3/8 LOCK WASHER |
| 19 | 4 | X-980783-03 | | | | #8 SST FLAT WASHER |
| 20 | 8 | X-980783-07 | | | | 3/8 SST FLAT WASHER |
| 21 | 28 | X-980794-02 | | | | 1/4 SST FLAT WASHER |
| 22 | 20 | X-980815-01 | | | | 1/4 INTERNAL TOOTH LOCK WASHER |
| 23 | 8 | X-980929-01 | | | | BOLT 3/8-16 X 2-3/4 SST |

| REV | ECO | DESCRIPTION | DATE | DFTM | CHKD |
|-----|-------|------------------|------------|------|------|
| E | 13580 | UPDATED HARDWARE | 12/28/2017 | TH | DSH |




* EXTENDED HEIGHT OPTION

BOLT THRU THESE (2) HOLES FOR 36" WORK SURFACE HEIGHT

BOLT THRU THESE (2) HOLES FOR 30" WORK SURFACE HEIGHT

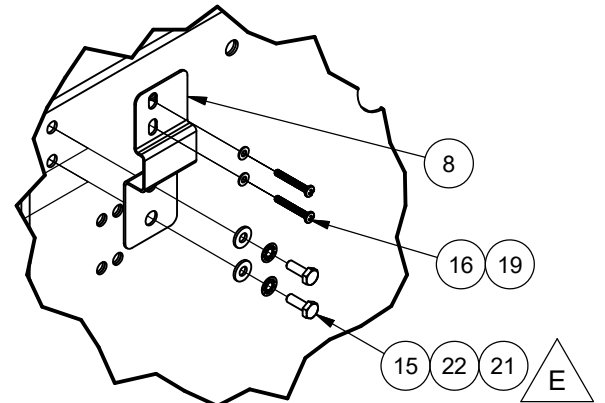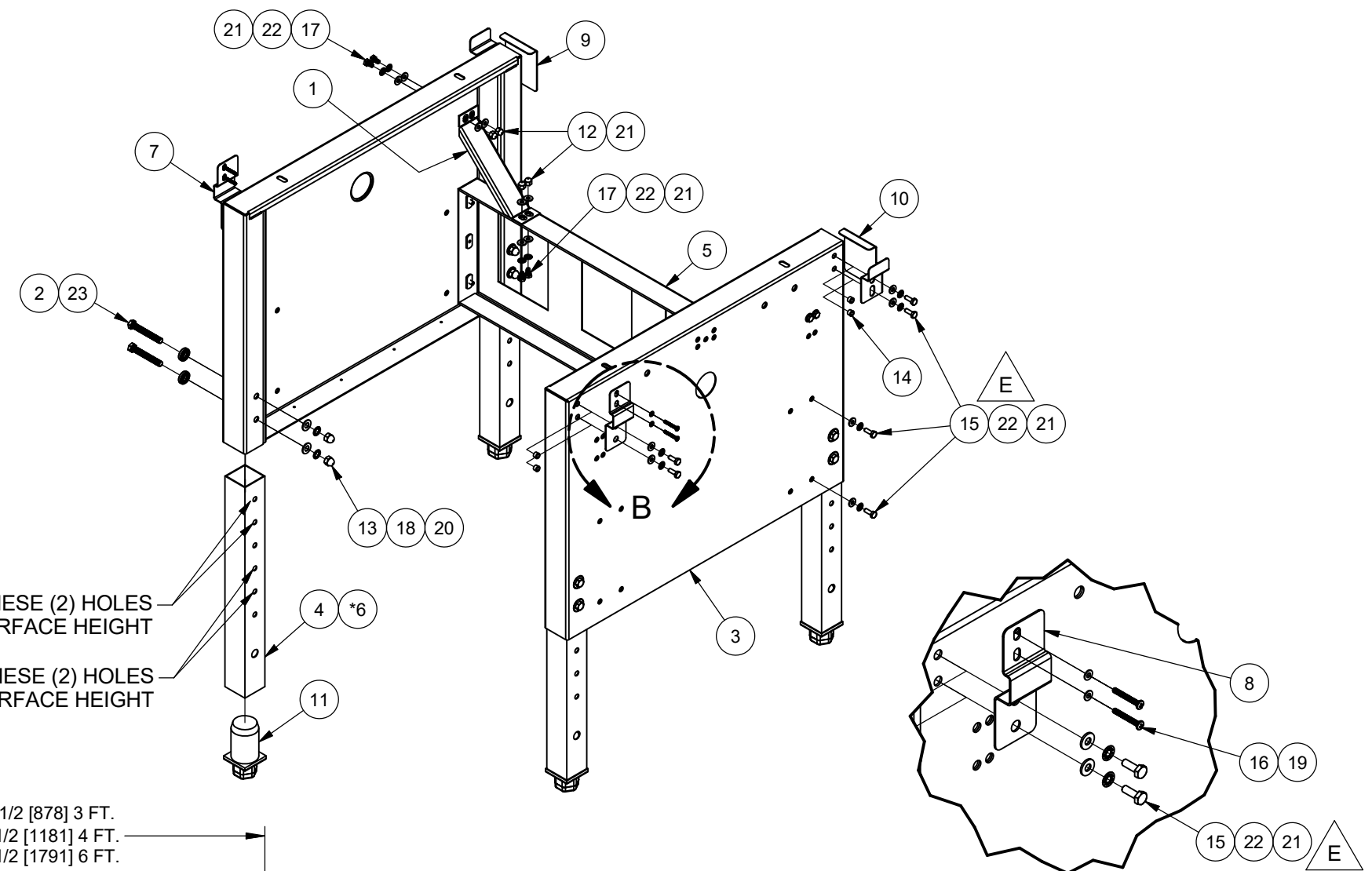

DETAIL B

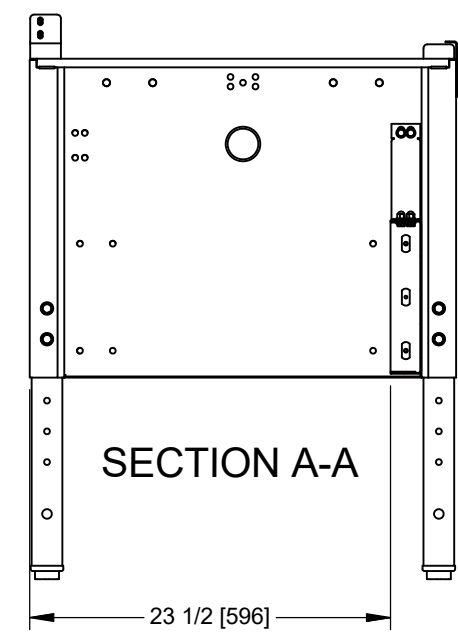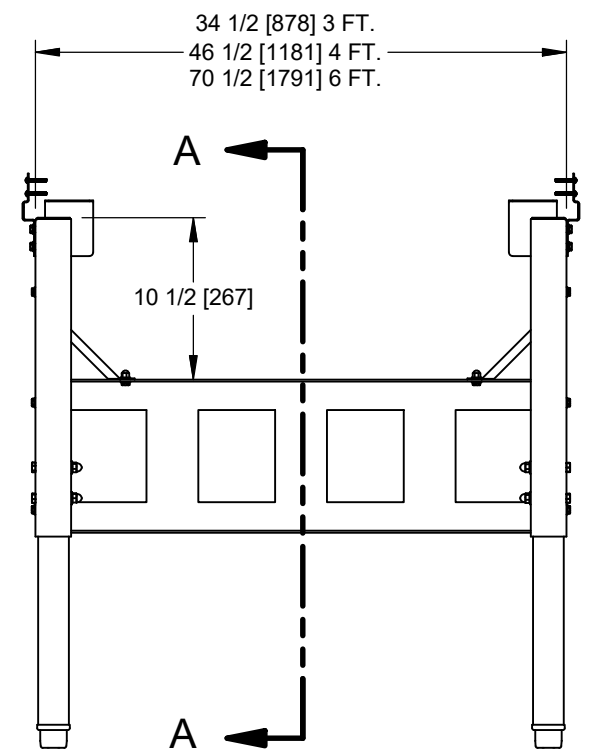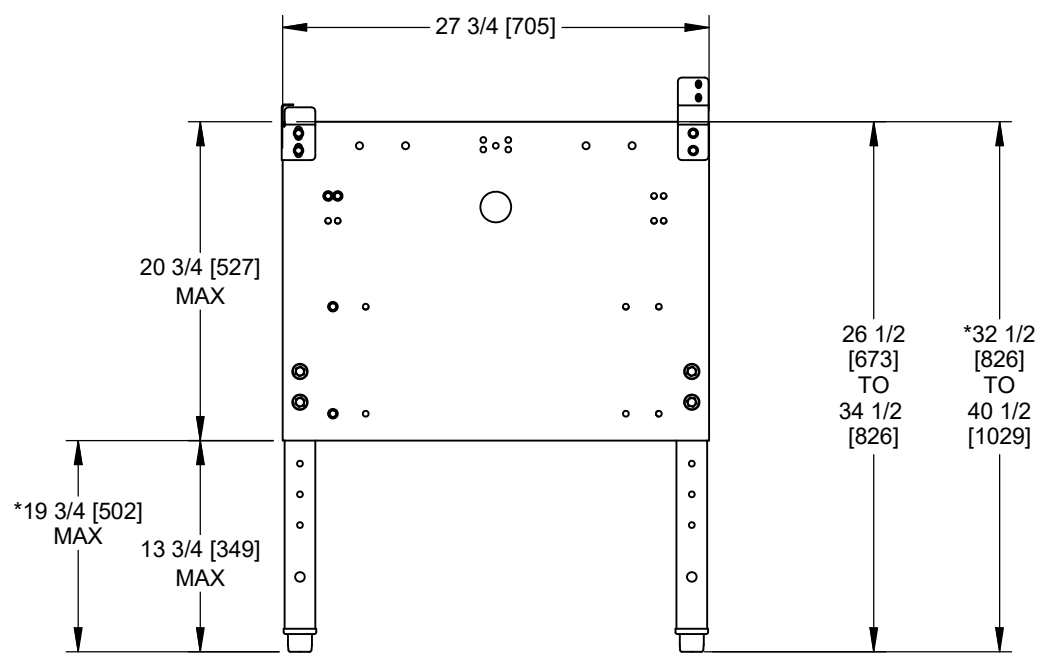

SECTION A-A

| E |
|------------|
| NU-813 |
| NU-800-128 |
| NU-800-129 |
| NU-800-130 |
| NU-800-140 |
| NU-800-141 |
| NU-800-142 |

| | | | | | |
|--|----------------|----------|-----------|------|---|
| | | | | | |
| TITLE ASSEMBLY, 28" BASESTAND TELESCOPING W/LL | | | | | |
| DFTM IW | DATE 4/13/2015 | CHKD MSS | SHEET 1 | OF 1 | |
| DRAWING NUMBER | | | BCD-17400 | | E |