

QUICK START GUIDE

Centrifugation Menu

Change Program Name

← N°10 → Program

← 5 → 7 → 15000 → 11:21 → 11

16000 RPM 20:00 min:sec 10°C

← Back to Home Screen

Press 5 Sec. to Lock or Unlock Program

← Acceleration / Braking Slopes

← Set Acceleration Slope

← Set Braking Slope

← Set Speed

← Set Timer

← Set Temperature (Refrigerated Model)

← Save Program

← Select Rotor

← Service Menu (Restricted Access)

← User Preferences

← Exit Back to Centrifugation Menu

← Hardware & Software Info.

← RPM RCF

← Greasing Reminder (in cycles)

← Show Usage Log

← Temp. Display Celsius or Fahrenheit

← Count At or Count Go

← Show Error Log

← Change Background Color

← End of Cycle Alarm ON

← End of Cycle Alarm OFF

← Run Refrigeration between Cycles

← Smart Pre-Cool

← Temp. Display Air or Sample

← Acceleration / Braking Slopes Graph

← Service Menu (Restricted Access) Details

← User Preferences Details

← Exit Back to Centrifugation Menu

← Hardware & Software Info.

← RPM RCF

← Greasing Reminder (in cycles)

← Show Usage Log

← Temp. Display Celsius or Fahrenheit

← Count At or Count Go

← Show Error Log

← Change Background Color

← End of Cycle Alarm ON

← End of Cycle Alarm OFF

← Run Refrigeration between Cycles

← Smart Pre-Cool

← Temp. Display Air or Sample

Home Screen

← Service Menu (Restricted Access)

← User Preferences

← Exit Back to Centrifugation Menu

← Hardware & Software Info.

← RPM RCF

← Greasing Reminder (in cycles)

← Show Usage Log

← Temp. Display Celsius or Fahrenheit

← Count At or Count Go

← Show Error Log

← Change Background Color

← End of Cycle Alarm ON

← End of Cycle Alarm OFF

← Run Refrigeration between Cycles

← Smart Pre-Cool

← Temp. Display Air or Sample

ON: timer starts when set speed is reached

OFF: timer starts when start is pressed